IKA

designed for scientists

IKA DRY BLOCK HEATER 1 IKA DRY BLOCK HEATER 2 IKA DRY BLOCK HEATER 3 IKA DRY BLOCK HEATER 4


Bezpečnostní pokyny	CS	48
Biztonsági utasítások	HU	49
Varnostna navodila	SL	50
Bezpečnostné pokyny	SK	51
Ohutusjuhised	ET	52
Drošības norādījumi	LV	53
Saugos nurodymai	LT	54
Указания за безопасност	BG	55
Instrucțiuni de siguranță	RO	56

Betriebsanleitung	DE	3
<small>Ursprungssprache</small>		
Operating instructions	EN	10
Mode d'emploi	FR	17
Инструкция по эксплуатации	RU	24
使用说明	ZH	31

Indicaciones de seguridad	ES	38
Veiligheidsaanwijzingen	NL	39
Norme di sicurezza	IT	40
Säkerhetsanvisningar	SV	41
Sikkerhedshenvisninger	DA	42
Sikkerhetsveiledning	NO	43
Turvallisuusohjeet	FI	44
Indicações de segurança	PT	45
Υποδείξεις ασφάλειας	EL	46
Wskazówki bezpieczeństwa	PL	47

**IKA DRY BLOCK HEATER 1
IKA DRY BLOCK HEATER 2
IKA DRY BLOCK HEATER 3
IKA DRY BLOCK HEATER 4**

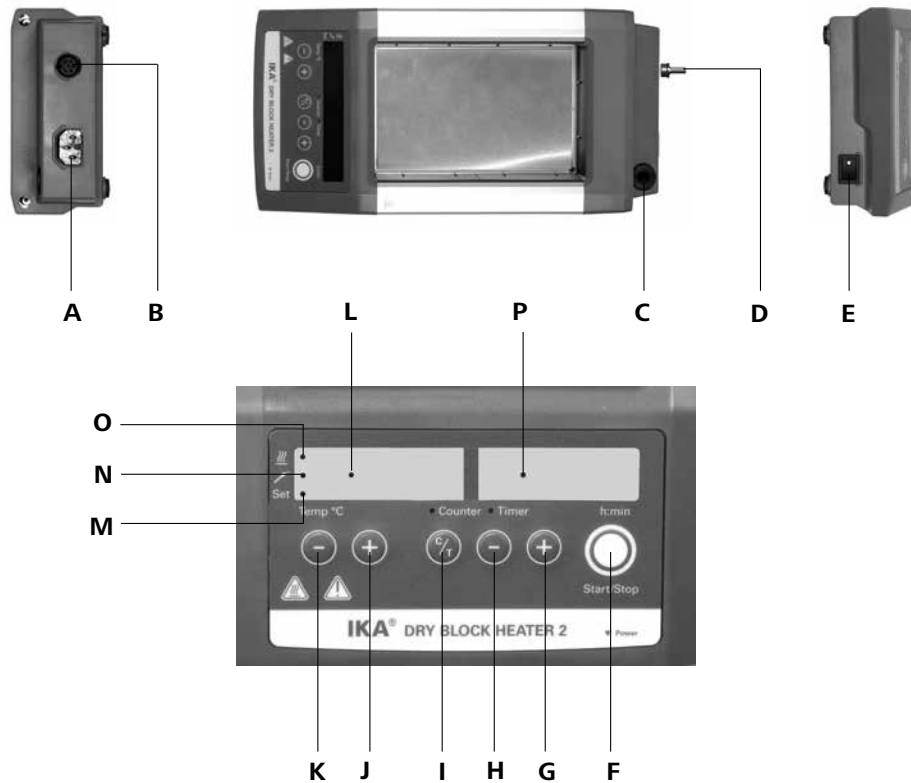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A	Netzbuchse
B	KT-Buchse
C	Stativgewindebohrung
D	Kontaktstecker
E	Geräteschalter
F	Taste "Start/Stop"
G	Taste Timer "+"
H	Taste Timer "-"
I	Taste Timer/ Counter
J	Taste Heizung "+"
K	Taste Heizung "-"
L	LED Display Heizung
M	LED (Set=Sollwert)
N	LED externer Temperatursensor
O	LED Heizplatte
P	LED Display Timer

Item Designation

A	Power socket
B	KT socket
C	Threaded hole for stand
D	Contact plug
E	Main switch
F	"Start/Stop" button
G	Button timer "+"
H	Button timer "-"
I	Button Timer/ Counter
J	Button heating "+"
K	Button heating "-"
L	LED Display heating
M	LED (set=set point value)
N	LED external temperature sensor
O	LED heating plate
P	LED display timer

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Contents

	Page
EU Declaration of conformity	10
Explication of warning symbols	10
Safety instructions	11
Intended use	12
Unpacking	12
Assembling the support rod	12
Operation	13
Setting the operating mode	13
Controlling the temperature via contact thermometer	14
Maintenance and cleaning	14
Error codes	15
Technical data	16
Warranty	16

EU Declaration of conformity

We declare under our sole responsibility that this product corresponds to the directives 2014/35/EU, 2014/30/EU and 2011/65/EU and conforms with the following standards or normative documents: EN 61010-1, EN 61010-2-010, EN 61326-1, EN 60529 and EN ISO 12100.

A copy of the complete EU Declaration of Conformity can be requested at sales@ika.com.

Explication of warning symbols

Indicates an (extremely) hazardous situation, which, if not avoided, will result in death, serious injury.

Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.

Indicates a potentially hazardous situation, which, if not avoided, can result in injury.

Indicates the exposure to a hot surface.

Safety instructions

General information

- Read the operating instructions in its entirety before using the device and follow the safety instructions.
- Keep the operating instructions in a place where it can be accessed by everyone.
- Ensure that only trained staff work with the device.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Electrical outlet must be grounded (protective ground contact).

Risk of burns! Exercise caution when touching parts of the housing and the heating plate.

The heating plate can reach dangerous temperatures. Pay attention to the residual heat on the heating plate after switching off the stirrer.

The device may only be transported when the heating plate has cooled down.

Device design

Do not use the device in explosive atmospheres, it is not EX-protected.

With substances capable of forming an explosive mixture, appropriate safety measures must be applied, e.g. working under a fume hood.

To avoid body injury and property damage, observe the relevant safety and accident prevention measures when processing hazardous materials.

- Set up the device in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- The feet of the device must be clean and undamaged.
- Ensure that the power cord set / temperature sensor cable does not touch the heating plate.
- Check the device and accessories for damage before each use. Do not use damaged components.

Permissible medium / contaminants / side reactions

Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.

- Beware of hazards due to:
 - flammable materials,
 - combustible media with a low boiling temperature,
 - glass breakage,
 - incorrect container size,
 - overfilling of media,
 - unsafe condition of container.
- Process pathogenic materials only in closed vessels under a suitable fume hood.

In unsupervised and safe operation, this device may only handle or heat media whose flash point is above the set safety temperature limit. (acc. to EN 61010-2-010)

- Please consider any possible contaminations and unwanted chemical reactions.

Procedures during sample runs

Wear your personal protective equipment in accordance with the hazard category of the media to be processed. There may

be a risk from:

- splashing and evaporation of liquids,
 - ejection of parts,
 - release of toxic or combustible gases.
- Do not heat liquid media except with the block. Never pour liquid media directly on to the heating plate!

Accessories

- Safe operation is guaranteed only with the use of original IKA accessories.
- Ensure that the external temperature sensor is inserted into the medium to a depth of at least 20 mm when connected.
- Always disconnect the plug before attaching accessories.
- Accessories must be securely attached to the device and cannot come off by themselves. The centre of gravity of the assembly must lie within the surface on which it is set up.
- Observe the operating instructions of the accessories.

Power supply / Switching off the device


The device will automatically restart in mode b following any interruption to the power supply.

- The voltage stated on the type plate must correspond to the mains voltage.
- The outlet for the mains plug must be easily accessible.
- The device can only be disconnected from the mains outlet by pulling out the power cord set or the connector plug.

For protection of the equipment

- The device may only be opened by qualified and IKA approved experts.
- Do not cover the device, even partially e.g. with metallic plates or film. This may result in overheating.
- Protect the device and accessories from bumps and impacts.
- Keep the base plate clean.
- Observe the minimum distances:
 - between devices min. 100 mm,
 - between device and wall min. 100 mm,
 - above the device min. 800 mm.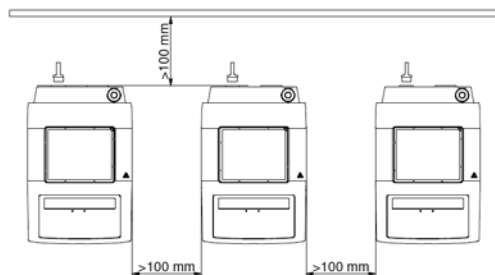

Intended use

• Use

- For heating media in block heaters

• Area of use

Indoor environments similar to that a laboratory of research, teaching, trade or industry area.

The safety of the user cannot be guaranteed:

- if the device is operated with accessories that are not supplied or recommended by the manufacturer,
- if the device is operated improperly or contrary to the manufacture's specifications,
- if the device or the printed circuit board are modified by third parties.

Unpacking

• Unpacking



- Please unpack the device carefully
- In the case of any damage a fact report must be sent immediately (post, rail or forwarder)

• Scope of delivery


- IKA DRY BLOCK HEATER
- Depending on the IKA DRY BLOCK HEATER model 1, 2, 3, or 4
- Mains cable
- Temperature sensor PT 1000.60
- Operating instructions

Assembling the support rod

- Review the mounting and safety instructions of the IKA boss head clamp prior to using it.
- The device must not be suspended from the support rod!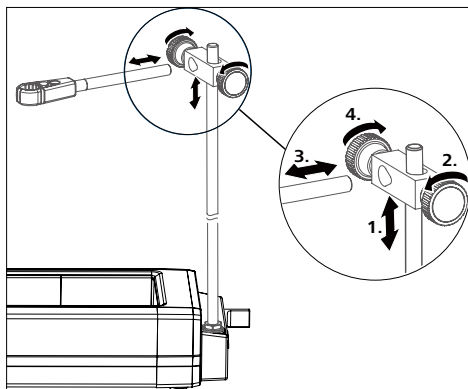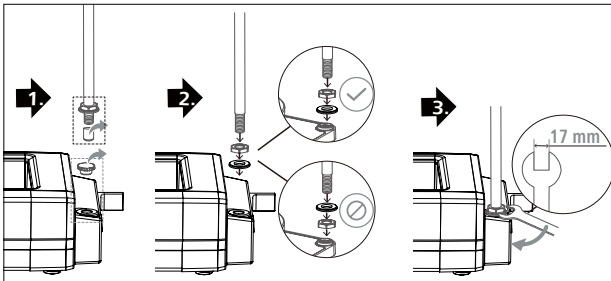

Risk of tipping! Make sure that the center of mass of the attached device does not protrude over the safe area indicated below with a dotted rectangle.

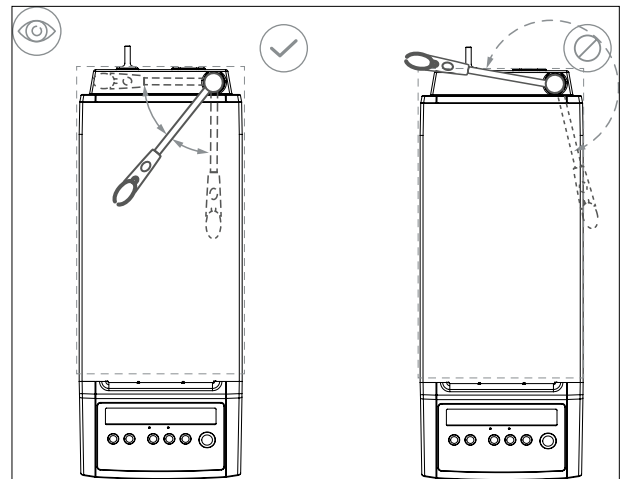Operation

	IKA DRY BLOCK HEATER 1	IKA DRY BLOCK HEATER 2	IKA DRY BLOCK HEATER 3	IKA DRY BLOCK HEATER 4
Commissioning	<ul style="list-style-type: none"> ☞ Move device switch (E) to the OFF position ☞ Insert the mains power cable into the power socket (A) ☞ Move device switch (E) to the ON position ⇒ The unit will be set to factory setting mode A (see "Setting operating mode") 			
Heating	<ul style="list-style-type: none"> ☞ Set the target temperature using the heater buttons +/- (J) and (K) ⇒ The value selected will be shown on the display (L) ☞ Start the heating function by pressing the "Start / Stop" button (F) 			
	<div style="border: 1px solid black; border-radius: 50%; width: 20px; height: 20px; display: flex; align-items: center; justify-content: center; margin: 5px;">i</div> <div style="border: 1px solid black; border-radius: 50%; width: 20px; height: 20px; display: flex; align-items: center; justify-content: center; margin: 5px;">i</div>	<p>The set-point and actual temperatures will be shown alternately on the display (L) is blinking: ⇒ When the heating is on, the LED (O) lights up.</p> <p>⇒ During standby operation, the display (L) will show HOT if the base plate temperature is above 50 °C</p>		
Connecting external thermometers (direct temperature control in the media)	<ul style="list-style-type: none"> ☞ Move device switch (E) to the OFF position ☞ Detach contact plug (D) ☞ Attach a DIN 12878 (Class 2) compliant safety contact thermometer or a PT 1000.60 temperature sensor to the socket (B) ☞ Move device switch (E) to the ON position 			
	<div style="border: 1px solid black; border-radius: 50%; width: 20px; height: 20px; display: flex; align-items: center; justify-content: center; margin: 5px;">i</div>	<p>Temperature sensor PT 1000 ⇒ The actual temperature for the temperature sensor shown on display (L) will correspond to the temperature of the media. The LED "external temperature sensor" PT 1000 (N) will be lit.</p>		

Setting operating mode

Operating the device in mode A or b

Mode A

All settings will be stored if the device is switched off or disconnected from the power supply. The heating functions will be set to OFF when the device is powered on.

Mode b

All settings will be stored if the device is switched off or disconnected from the power supply. The heating function will be set to ON or OFF when the device is powered on, depending on the previous status of the device.

Factory setting: mode A

The mode selected will be shown on the display when the device is started up.

Change the mode

- ☞ Move device switch **(E)** to the OFF position
- ☞ Keep the selection button **(F)** depressed
- ☞ Move device switch **(E)** to the ON position
- ☞ Release the selection button **(F)**

i

 ⇒ The set mode is indicated on the Display **(P)** b

Timer / Counter

After the device has been switched on, the "Timer/ Counter" button **(I)** can be used to toggle between timer mode and counter mode.

Timer

Pressing the buttons "+" **(G)** and "-" **(H)** allows selection of a time period of up to 99 hours and 59 minutes. Once the heating function **(F)** is activated, the preset time starts to run. When the total time has run down to 00 hours and 00 minutes, a warning buzzer sounds for 30 seconds and in addition the display flashes. The heating function is then deactivated.

Counter

When the heating function is stated by pressing the button **(F)** the counter is activated. In counter mode the time is counted upwards from 00 hours 00 minutes until the heating function is stopped. The counter always starts from 00:00 hh:mm.

Selection of Timer / Counter

Pressing the button **(I)** toggles between timer mode and counter mode. The mode can be activated only when the heating function is off.

Controlling the temperature via contact thermometer

The preferable method for controlling the average temperature is with the contact thermometer. After the set point temperature has been adjusted, this results in a short heating-up time, practically no temperature drift and only minor fluctuation in temperature.

A 6-pin jack is located on the rear side of the instrument for connecting the contact thermometer or the contact plug. The electronics of the instruments returns a test current that must flow via connector pins 3 and 5 for the heating plate to heat up.

Settings

For the detailed instructions for settings and limit values, refer to the operating instructions for the instrument you are connecting.

The target temperature of the medium is set at the contact thermometer. The heater buttons +/- **(J) and **(K)** on the device can be used to preselect the required surface temperature of the heating plate.**

Safety contact thermometers

acc. to DIN 12 878 class 2 or acc. Gerstel are connected with a three-wire cable, the test current flows through the contact thermometer.

Safety function

If the test current is interrupted because of e.g. breakage of contact thermometer or falling out of the cable plug, the heating cuts off. The required plug allocation is indicated on a reference plate above the diode bushing.

Contact thermometer without safety circuit

acc. to DIN 12 878 class 0. The instrument only heats if the test current circuit is closed by an electrical connection of the plug pins 3 and 5.

Maintenance and cleaning

The device is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

Cleaning

- For cleaning disconnect the mains plug!
- Use only cleaning agents which have been approved by IKA to clean IKA devices.
These are water (with tenside) and isopropanol.
- Wear protective gloves during cleaning the devices.
- Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.
- Do not allow moisture to get into the device when cleaning.
- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.

Spare parts order

When ordering spare parts, please give:

- device type.
- serial number, see type plate.
- position number and description of spare part, see **www.ika.com**.
- software version

Repair

Please only send devices in for repair that have been cleaned and are free of materials which might present health hazards.

For repair, please request the "Safety Declaration (Decontamination Certificate)" from IKA or use the download printout of it from IKA website at **www.ika.com**.

If your appliance requires repair, return it in its original packaging. Storage packaging is not sufficient when sending the device - also use appropriate transport packaging.

Error codes

Any malfunctions during operation will be identified by an error message on the display (LED **(P)**).

Proceed as follows in such cases:

- Switch off device using the main switch **(E)**
- Carry out corrective measures
- Restart device

Error code	Cause	Effect	Solution
E 3	Temperature inside device too high	Heating off	Switch off device and allow to cool down
E 6	Break in safety circuit	Heating off	<ul style="list-style-type: none"> - Plug in contact plug (D) - Plug in PT 1000 contact thermometer / temperature sensor - Replace faulty connecting cable, plug, or contact thermometer
E 25	Heating and switching element monitoring	Heating off	<ul style="list-style-type: none"> - Switch off device - Warning! Only to be carried out by authorised service personnel: Carry out an internal test on the device to check the plug-in connector for the heating element
E 26	Difference between temperature of safety sensor and temperature of control sensor control temperature > (safety temperature + 40 K)	Heating off	<ul style="list-style-type: none"> - Switch off device - Warning! Only to be carried out by authorised service personnel: Carry out an internal test on the device to check the temperature sensor

If the actions described fail to resolve the fault or another error code is displayed then take one of the following steps:

- contact the **IKA** service department,
- send the device for repair, including a short description of the fault.

Technical Data

		IKA DRY BLOCK HEATER 1	IKA DRY BLOCK HEATER 2	IKA DRY BLOCK HEATER 3	IKA DRY BLOCK HEATER 4
Operating voltage range	Vac	(220 ... 230) ± 10%			
	Vac	115 ± 10%			
	Vac	100 ± 10%			
Rated voltage	Vac	230 / 50 Hz			
	Vac	115 / 60 Hz			
	Vac	100 / 60 Hz			
Frequency	Hz	50 / 60			
Display		digital			
Temperature range	°C	ambient temperature +5 ... 120			
Timer	min/h	1 min ... 99 h 59 min			
Perm. duration of operation	%	100			
Perm. ambient temperature	°C	+5 ... +40			
Perm. humidity	%	80			
Protection acc. to DIN EN 60529		IP 21			
Protection class		I			
Overvoltage category		II			
Contamination level		2			
Operation at a terrestrial altitude	m	max. 2000			
Dimension (W x D x H)	mm	151 x 228 x 73	151 x 304 x 73	151 x 380 x 73	151 x 456 x 73
Weight	kg	1.3	1.7	2.0	2.0
Heating					
Power consumption (+10%) max. at 230 Vac	W	165	250	330	412
115 Vac		165	250	330	412
100 Vac		165	250	330	412
Temperature sensor PT 1000 variation DIN EN 60751 Kl. A	K	≤ ± 0.1*			
Temperature stability within the blocks at 37 °C	°C	± 0.2	± 0.2	± 0.2	± 0.2
Temperature stability within the blocks at 60 °C	°C	± 0.4	± 0.4	± 0.4	± 0.4
Temperature homogeneity within the the blocks at 37 °C	K	± 0.2	± 0.2	± 0.3	± 0.3
Temperature homogeneity within the the blocks at 60 °C	K	± 0.4	± 0.4	± 0.6	± 0.6
Heating rate (PT 1000 in solid block)	K/min	5.0	4.5	4.0	4.0
Fixed safety circuit					
Safety temperature limit	°C	150			

* Test parameter: with test block and glass vessel filled with M50 oil

Subject to technical changes!

Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited