2000004385

C 1.30_112015

C 1.30

Betriebsanleitung	DE	3
Ursprungssprache		
Operating instructions	EN	6
Mode d'emploi	FR	10
Руководство пользователя	RU	13

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Contents

EN

	Page		Page
Declaration of conformity	2	Function	8
Contents	6	Maintenance and cleaning	9
Explanations of symbols	6	Fault table	9
Safety instructions	6	Warranty	9
Unpacking	7	Spare part list	17
Setting up of the instrument	8	Spare parts	18

Explanations of symbols

General hazard

This symbol identifies information **which is of absolute importance for the safety of your health**. Non-compliance may lead to impairments to health and injury.

This symbol identifies information which is important for proper operation and handling of the decomposition vessel. Non-compliance may result in imprecise measurements or damage to the venting station.

Safety instructions

Read the operating instructions in full before starting up and follow the safety instructions.

Keep the operating instructions in a place where they can be accessed by everyone.

Ensure that only trained staff work with the appliance.
Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

Furthermore, toxic combustion residue in the form of gases, ashes or condensation, for example, is possible in the inner wall of the washing bottle.

⚠ DANGER Observe the accident prevention requirements applicable to the activity and the work station.

Wear your personal protective equipment.

When handling combustion samples, combustion residue and auxiliary materials, please observe the relevant safety regulations. The following materials, for example, could pose a risk:

- corrosive
- highly flammable
- explosive
- bacteriologically contaminated
- toxic

⚠ DANGER Combustion gases are hazardous to health, therefore the venting hose must be connected to a suitable gas cleaning system or extraction system.

⚠ DANGER The C 1.30 venting station is used for controlled venting of IKA C1 calorimeter.

The venting speed may be infinitely adjusted via a throttle valve by using a manual adjusting screw. Please note that the throttle valve is **closed** at the beginning of the venting operation.

Do not operate without the ground clamp included in the delivery quantity!

The maximum permitted operating pressure is 40 bar (at the pressing tubing **(d)**).

Please observe the general laboratory guidelines when handling glass vessels.

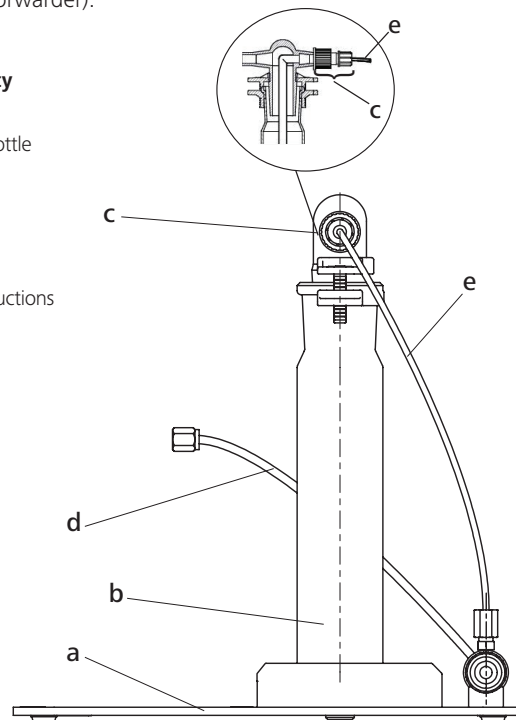
Unpacking

• Unpacking

- Please unpack the device carefully
- In the case of any damage a report must be sent immediately (post, rail or forwarder).

• Delivery quantity

- Base plate
- Gas washing bottle
- GL screw joint
- Pressure tubing
- Drainage hose
- Fork spanner
- Set of O-Rings
- Operating instructions

Setting up in the instrument

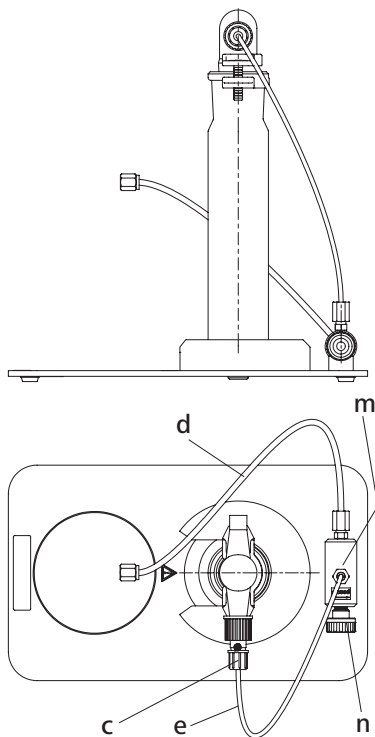
Place the venting station on a horizontal surface and insert the gas washing bottle into the gas washing bottle holder until it comes to a stop. Connect the drainage hose (e) to the throttle valve (d) and the GL screw joint (c) of the washing bottle as shown in the diagram. The tubing is fastened by closing the GL screw joint.

WARNING

Now fasten the pressure tubing (d) to the C 1 calorimeter. Dis-

connect the flexible venting tube from the screw SW 10. The capillary and the screw remain at C 1.

All of the screw joints are tightened by means of the SW 10 fork spanner (1/4 turning) which is included in the delivery quantity.

Function

Pour an absorption solution that is suited to your purposes into the gas washing bottle and **secure the ground joints by means of the ground clamp**. Close the throttle valve by slightly turning the adjusting screw in a clockwise direction.

Note:

There may be some formation bubbles in the system when the adjusting screw is hand-tightened with maximum operating pressure in the decomposition vessel.

Check to make certain that the ground clamp fits correctly onto the gas washing bottle and slowly open the throttle valve.

Take note of any bubble formation and correct them in order to achieve the venting speed desired.

If the C 1 is completely vented and the water is drained, it can be opened.

WARNING

The gas washing bottle may be damaged if venting is performed too quickly.

Set on the C 1 in the menu "Basic adjustment" the value for "Extend. O₂ Emptying" to 5 min. or longer. Make sure that the pressure in C 1 within the set time is emptied. Do this, turn the throttle screw (n) according to.

Maintenance and cleaning

Only use cleansing agents which have been recommended by **IKA**®.

Use to remove:

Dyes	Isopropyl alcohol
Construction materials	Water containing surfactant, Isopropyl alcohol
Cosmetics	Water containing surfactant, Isopropyl alcohol
Foodstuffs	Water containing surfactant
Fuels	Water containing surfactant

- Do not allow moisture to get into the appliance when cleaning.
- Wear protective gloves when cleaning the devices.
- Please consult with IKA before using any cleaning or decontamination methods, other than those recommended here.

Ordering spare parts

When ordering spare parts, please give:

- Device type
- Manufacturing number, see type plate
- Item number and designation of the spare part, see www.ika.com.

Repair

Please send instrument in for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.

For this you should request the **"Decontamination Clearance Certificate"** from **IKA**®, or use the download printout of it from the **IKA**® website www.ika.com.

Return the instrument in its original packaging. Storage packaging is not sufficient. Also, please use suitable shipping package materials.

Fault table

Error:	C 1 is not expanded.
Cause:	Capillary at C1 blocked. Throttle valve does not open.
Remedy:	Replace capillary. Open throttle screw.
<hr/>	
Error:	Leakage on throttle valve
Cause:	O-ring on throttle screw defective.
Remedy:	Replace O-ring.
<hr/>	

Warranty

In accordance with **IKA**® warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine directly to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

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IKA® - Werke GmbH & Co.KG