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EN

Useful information

Observe the operating instructions of the calorimeter C 1.

For the sample preparation use the Organizer C 1.1012 (delivery scope calorimeter C1).

The combustion chamber C 1.10 consists of stainless steel.

The halogen resistant combustion chamber C 1.12 (Hastelloy C 4) has been designed for anorganic samples with higher halogen content, particularly chlorine samples.

To enable the quantitative evaluation of halogens and sulphur, the combustion chamber C 1.12 is preconditioned before shipment. This means that the inner surfaces and the fittings of the combustion chamber are catalytically activated by means of a patented procedure. Parts thus treated are recognizable because of their opaque, tarnished and partly spotty surfaces. This appearance of this surface is intentional and should not be confused with traces of wear.

This state is important for the following analysis and may not be treated with brushes or other hard materials. Wiping and drying should be done using a soft cloth.

The combustion Chamber C 1.10 can not be used for such applications because of the safety risk and insufficient analysis results.

Safety instructions

Observe the operating instructions of the calorimeter C 1. Read the operating instructions completely before starting up and follow the safety instructions.

- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the equipment.

WARNING

Check the seals for damage before each use.

CAUTION

Observe the accident prevention requirements applicable to the activity and the work station. Wear your personal protective equipment. When handling combustion samples, combustion residue and auxiliary materials, please observe the relevant safety regulations.

DANGER

Do not incinerate metal-containing samples.

- The following materials, for example, could pose a risk: corrosive, highly flammable, explosive, bacteriologically contaminated, toxic.

Correct use

The combustion chambers C 1.10 and C 1.12 may only be operated with the calorimeter C 1.

The combustion chamber C 1.12 is recommended, if the samples have unknown ingredients or higher chlorine content (> 1 %), e.g. PVC or halogenated solvents.

Main application field: disposal industry

Unpacking

- **Unpacking**
 - Please unpack the device carefully
 - In the case of any damage a detailed report must be sent immediately (post, rail or forwarder).

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- Verbrennungskammer
- C 1.103 Zünddraht (C 1.10)
- C 1.123 Platin-Zünddraht (C 1.12)
- C 710.4 Baumwollfäden
- C 6 Quarztiegel
- Griff
- O-Ring (3,1 x 1,6)
- O-Ring (48 x 2,0)
- Sechskantmutter M3 Hastelloy
- Doppelmaulsschlüssel (5,5 x 7,0)
- Doppelmaulsschlüssel (8,0 x 10,0)

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The ignition wire must be fixed- if necessary tighten the hexagon nut (M3 Hastelloy) with the double open ended spanner.

After each combustion are all parts of the combustion chamber that come in contact with the combustion products to clean condensate. It is sufficient to parts with an absorbent, lint-free cloth to wipe off.

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	C 1.10	C 1.12
Werkstoff	Edelstahl	Hastelloy
Volumen	160 ml	160 ml
Zünddraht	Kanthal	Platin


Technical data

	C 1.10	C 1.12
Material	Sainless steel	Hastelloy
Volume	160 ml	160 ml
Inginition wire	Kanthal	Platinum

IKA®-Werke GmbH & Co.KG

Janke & Kunkel-Str. 10
 D-79219 Staufen
 Tel. +49 7633 831-0
 Fax +49 7633 831-98
 sales@ika.de

www.ika.com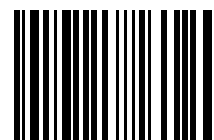

20011171