IKA

designed for scientists

IKA HELISTAND

5

/// ENGLISH

15

/// SVENSKA

19

/// NORSK

23

/// ČEŠTINA

27

/// SLOVENSKI

31

/// EESTI

35

/// LIETUVOS

39

/// ROMÂNĂ

13

/// NEDERLANDS

17

/// DANSK

21

/// SUOMI

25

/// MAGYAR

29

/// SLOVAK

33

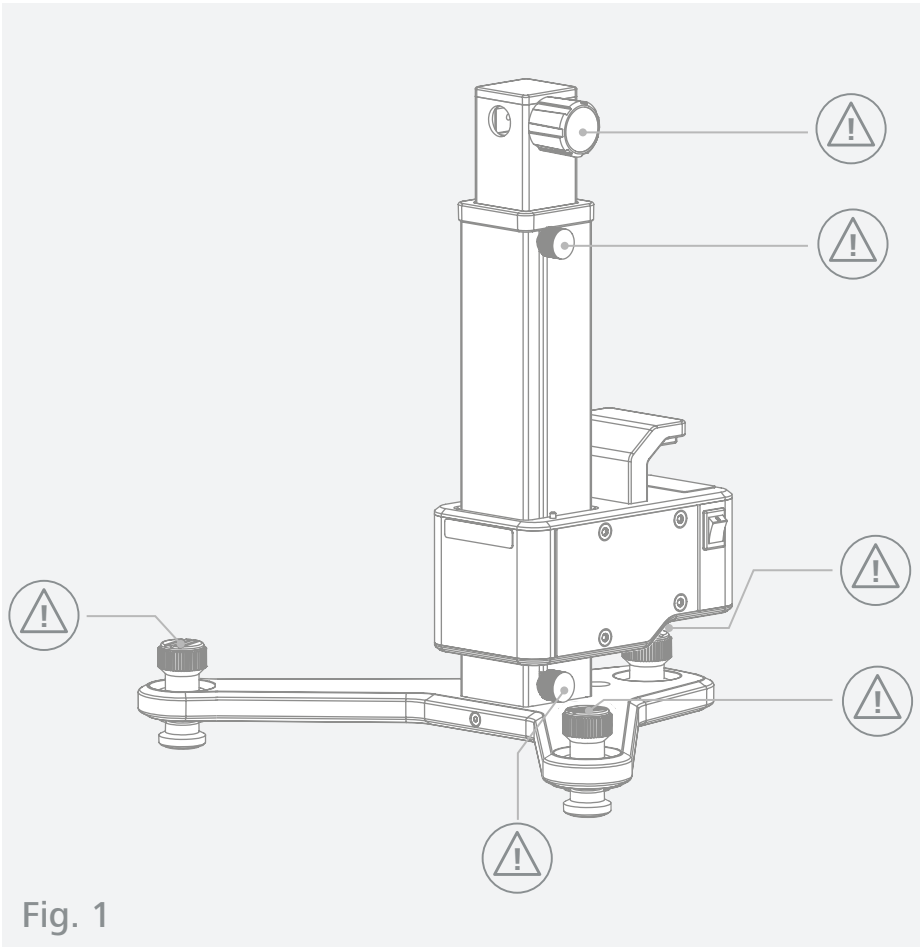
/// LATVISKI

37

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41

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	EU Declaration of conformity	5
	Explication of warning symbols	6
	Safety instructions.....	6
	Intended use.....	7
	Unpacking	8
	Installation.....	8
	Operation	9
	Maintenance and cleaning	12
	Technical Data	12
	Warranty	12

Source language: German

EU Declaration of conformity

We declare under our sole responsibility that this product corresponds to the directives 2014/35/EU, 2006/42/EC, 2014/30/EU and 2011/65/EU and conforms with the following standards or normative documents: EN 61010-1, EN 61326-1, EN 60529 and EN ISO 12100. A copy of the complete EU Declaration of Conformity can be requested at sales@ika.com.

Explication of warning symbols

/// Warning symbols

Caution! Indicates a potentially hazardous situation, which, if not avoided, can result in injury.

Notice! Indicates practices which, if not avoided, can result in equipment damage.

/// General Symbols

A — Position number
Indicates device components relevant to actions.

Correct / result
Indicates the correct execution or the result of an action step.

Wrong
Indicates the incorrect execution of an action step.

Note
Indicates steps of actions that require particular attention.

Safety instructions

/// General information

- › **Read the operating instructions in its entirety before using the device and follow the safety instructions. If this accessory is used with another device, observe also its operating instructions.**
- › Keep the operating instructions in a place where it can be accessed by everyone.
- › Ensure that only trained staff work with the device.
- › Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- › Electrical outlet must be grounded (protective ground contact).

Notice!

- › Pay attention to the marked sites in Fig. 1.

/// Device design

Notice!

- › Set up the device in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- › The feet of the device must be clean and undamaged.
- › All screw connections must be properly tightened.
- › Check the mounting position of the viscometer to avoid the danger of the unit tipping over.

/// Working with the device

Notice!

- › When transporting the stand, ensure the device is in its lowest (end stop) position.

/// Accessories

- › Protect the device and accessories from bumping and impacting.
- › Check the device and accessories beforehand for damage each time when you use them. Do not use damaged components.
- › Safe operation is guaranteed only with the use of original IKA accessories.
- › Always disconnect the plug before attaching accessories.

/// Power supply / Switching off the device

- › The voltage stated on the type plate must correspond to the mains voltage.
- › The outlet for the mains plug must be easily accessible.
- › The device can only be disconnected from the power supply by pulling out the power plug or the connector plug.

Notice!

- › When lifting and lowering, make sure that the power cable can move freely.

Intended use

/// Use

- › The IKA HELISTAND was designed for comparative measurements for samples like creams, pastes, gels by lifting and lowering and for holding the IKA ROTAVISC. The automatic up and down movement moves the T-spindle mounted on the IKA ROTAVISC helically through the sample, therefore the channeling problem will be eliminated.

Notice!

- Operation at low speeds will usually produce optimum readings. Recommended speed:
 - lo-vi: max. 12 rpm
 - me-vi, hi-vi I and hi-vi II: max. 10 rpm
- Activate the IKA HELISTAND drive only when using the T-spindle.
- The IKA HELISTAND can only be used without the IKA ROTAVISC guard rail.
- The IKA HELISTAND (motorized mode) is not recommended for checking the calibration of the IKA ROTAVISC.

/// Area of use

- › Indoor environments similar to that a laboratory of research, teaching, trade or industry area.
- › The safety of the user cannot be guaranteed:
 - if the device is operated with accessories that are not supplied or recommended by the manufacturer,
 - if the device is operated improperly or contrary to the manufacture's specifications,
 - if the device or the printed circuit board are modified by third parties.

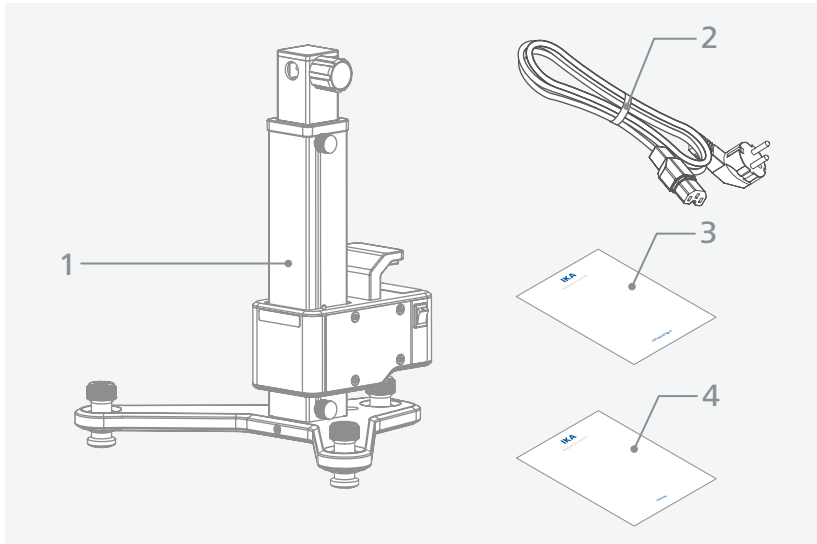
Unpacking

/// Unpacking

- › Unpack the device carefully. Any damage should immediately be reported to the carrier (mail, rail or freight forwarding company).

/// Scope of delivery

1	IKA HELISTAND
2	Power cord
3	Operating instructions
4	Warranty card

Installation

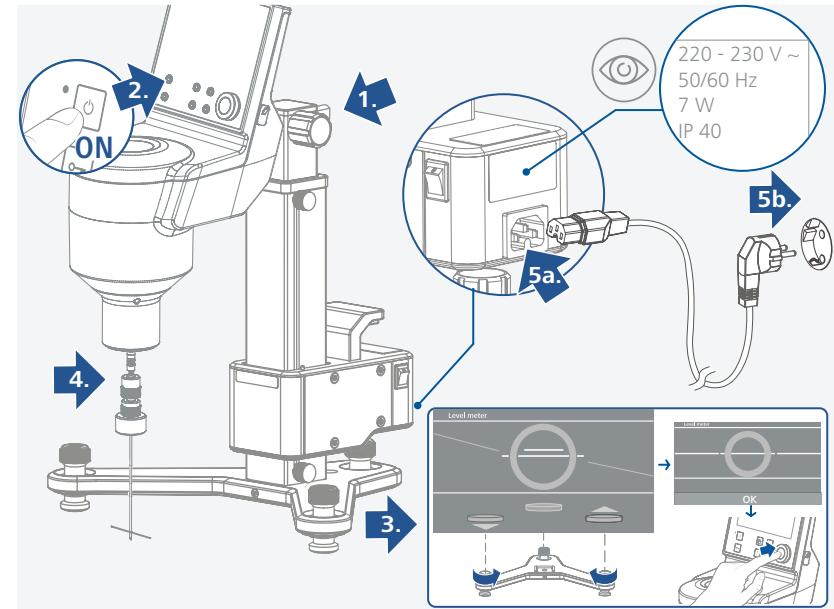
/// Assemble IKA ROTAVISC to IKA HELISTAND

- › Refer to the operating instructions of IKA ROTAVISC.

/// Assemble the spindle

- › Refer to the operating instructions of T-SP-Set.


/// Connect IKA HELISTAND to a socket


Operation

/// Height adjustment

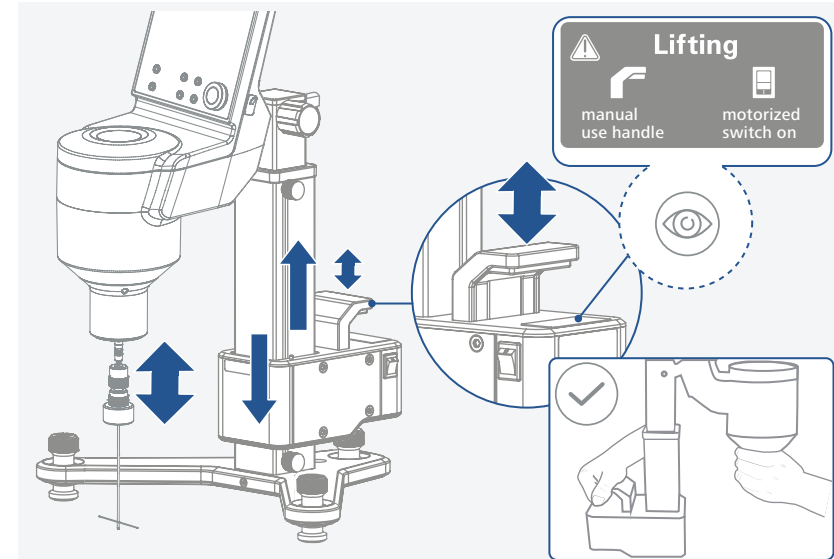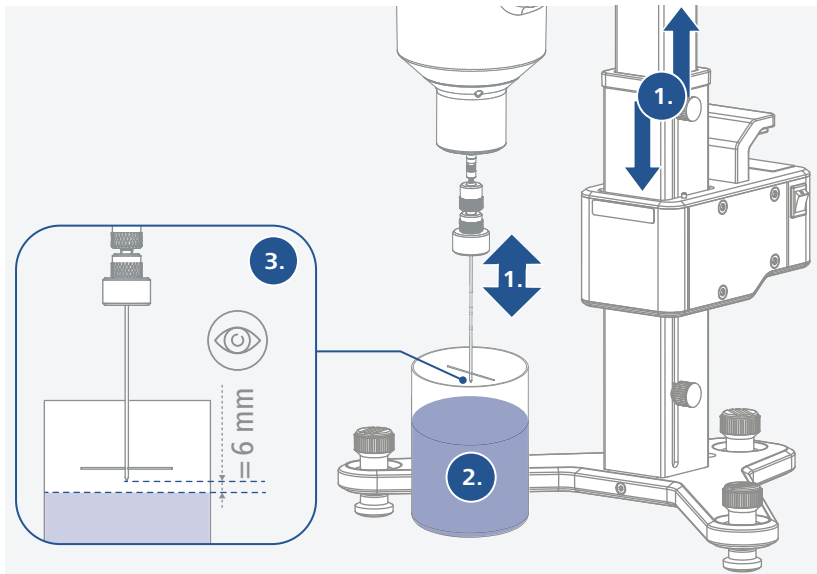
- › Adjust the spindle initial position while unlocking and lifting the handle.
- › The initial position of the spindle peak should be about 6 mm above the sample level.


Caution!

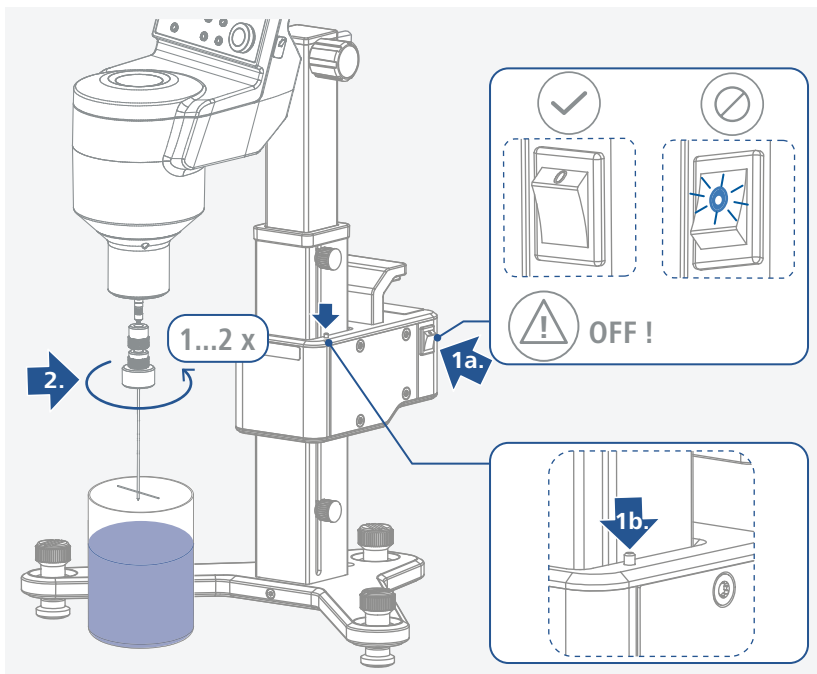- › Make sure that the IKA HELISTAND is switched off!
- Note:** The switch indicator light is off.
- › Push the pin downwards, the IKA HELISTAND will move down when the drive is switched on.
- › Let the spindle run for one to two revolutions before turning on the IKA HELISTAND.


/// Set limit stops

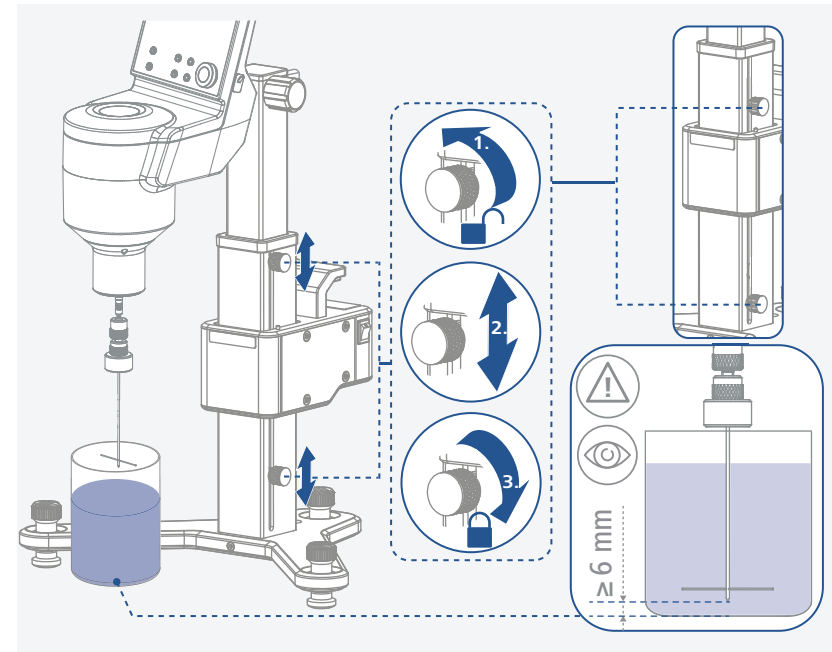- › Adjust the lower and upper limit stops. Make sure that the adjustment of the limit stops corresponds to the required penetration depth of the spindle.
- Note:** The minimum allowed distance between spindle and vessel bottom is 6 mm.

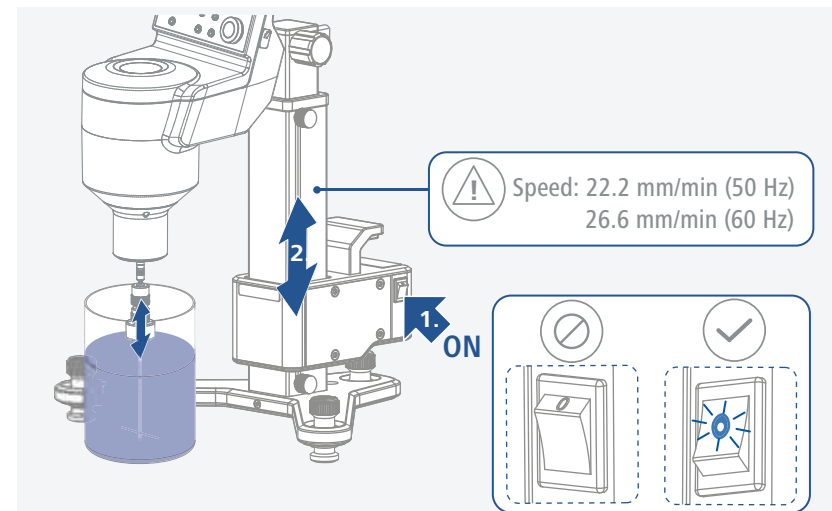- › Switching on: Switch on the drive by flipping the switch.
- Note:** The switch indicator light will illuminate.
- The speed of the up-and-down movement is 22.2 mm/min.

Maintenance and cleaning

- › The device is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

/// Cleaning

- › For cleaning disconnect the mains plug!
- › Use only cleaning agents which have been approved by IKA to clean IKA devices. These are water (with tenside) and isopropanol.
- › Wear protective gloves during cleaning the devices.
- › Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.
- › Do not allow moisture to get into the device when cleaning.
- › Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.

/// Spare parts order

- › When ordering spare parts, please give:
 - device type.
 - serial number, see type plate.
 - position number and description of spare part, see www.ika.com.

/// Repair

- › Please only send devices in for repair that have been cleaned and are free of materials which might present health hazards.
- › For repair, please request the "Safety Declaration (Decontamination Certificate)" from IKA or use the download printout of it from IKA website at www.ika.com.
- › If your appliance requires repair, return it in its original packaging. Storage packaging is not sufficient when sending the device - also use appropriate transport packaging.

Technical data

Voltage	220 ... 230 VAC ± 10%
	100 ... 115 VAC ± 10%
Frequency	50 / 60 Hz
Power input	7 W
Power input max.	7 W
Permissible ambient temperature	+ 5 ... + 40 °C
Permissible relative humidity	80 %
Protection class according to DIN EN 60529	IP 40
Overload protection	yes
Protection class	I
Contamination level	2
Overvoltage category	II
Operation at a terrestrial altitude	max. 2000 m above sea level
Stroke max.	142 mm
Diameter for extension rod (Ø)	16 mm
Dynamic load	3,2 kg
Height adjustable safety switch	yes
Dimensions (W x D x H)	351 x 252 x 343 mm
Weight	5,5 kg

Subject to technical changes!

Warranty

- › In accordance with IKA warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.
- › The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited