IKA

designed for scientists

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Contents

EN

| | Page |
|--|------|
| EU Declaration of conformity | 11 |
| Safety instructions | 11 |
| Intended use | 12 |
| Unpacking | 12 |
| Commissioning | 13 |
| Starting | 13 |
| Function stirring | 13 |
| Function heating | 14 |
| Limiter of safety temperature | 14 |
| Adjusting the Safety Temperature limit "Safe Temp" (Vers. KT) | 15 |
| Controlling the Medium Temperature via ETC 1 or Contact Thermometer (Vers. KT) | 15 |
| Error codes | 16 |
| Mounting - Support H 16 V | 16 |
| Maintenance and cleaning | 17 |
| Accessories | 17 |
| Warranty | 17 |
| Technical data | 18 |

EU Declaration of conformity

We declare under our sole responsibility that this product corresponds to the directives 2014/35/EU, 2006/42/EC, 2014/30/EU and 2011/65/EU and conforms with the following standards or normative documents: EN 61010-1, EN 61010-2-010, EN 61010-2-051, EN 61326-1, EN 60529 and EN ISO 12100.

A copy of the complete EU Declaration of Conformity can be requested at sales@ika.com.

Safety instructions

The **IKA RH basic 2; RH basic KT/C or RH digital KT/C** was not constructed for operations in dangerous atmospheres, for mixing dangerous substances or for operation under water.

- › Do not use the device in explosive atmospheres, it is not EX-protected.
- › With substances capable of forming an explosive mixture, appropriate safety measures must be applied, e.g. working under a fume hood.
- › To avoid body injury and property damage, observe the relevant safety and accident prevention measures when processing hazardous materials.
- › Set up the device in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.

Attention – Magnetism! Effects of the magnetic field have to be taken into account (e.g. data storage media, cardiac pacemakers ...).

- › Ensure that the power cord set / temperature sensor cable does not touch the heating plate. The power cable must be replaced by a cable of equal quality and capacity.
- › **Version RH basic KT/C and RH digital KT/C:** Examine regularly (we recommend every 4 weeks) the function of the safety temperature limiter. For this the safety temperature must be changed around at least 1/4 revolution of the value already stopped and be accomplished the attitude as under "adjusting the safety temperature delimitation SAFE temp" described. By the exceeding of the adjusted safety temperature (e.g. rotary button "temp" posed on right end stop) you can examine the function. The "temp" pilot lamp must expire and the heating must be lasting switched off.
- › When using PTFE-coated magnetic bars, the following has to be noted: Chemical reactions of PTFE occur in contact with molten or solute alkali metals and alkaline earth metals, as well as with fine powders of metals in groups 2 and 3 of the periodic system at temperatures above 300 °C – 400 °C. Only elementary fluorine, chlorotrifluoride and alkali metals attack it; halogenated hydrocarbons have a reversible swelling effect.

(Source: Römpps Chemie-Lexikon and "Ulmann", Volume 19)

Risk of burns! Exercise caution when touching parts of the housing and the heating plate. The heating plate can reach dangerous temperatures. Pay attention to the residual heat on the heating plate after switching off the stirrer.

The device may only be transported when the heating plate has cooled down.

- › Please consider any possible contaminations and unwanted chemical reactions.
- › Before placing the device in service, please ensure that the turn dial for adjusting the speed is against the left stop, since the device will begin to run at the last speed that was set.
- › The device will automatically restart following any interruption to the power supply.
- › The device may not be transported by using an H 16 V support rod or the power cable.
- › Clicking noises may occur as a result of material expansion during heating and cooling.
- › In unsupervised and safe operation, this device may only handle or heat media whose flash point is above the set safety temperature limit. (acc. to EN 61010-2-010)
- › Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- › The device may only be opened by qualified and IKA approved experts.
- › Safe operation is guaranteed only with the use of original IKA accessories.

Intended use

Use

- › The IKA RH basic 2; RH basic KT/C or the IKA RH digital KT/C is a stirring instrument with heating function. The instrument is suitable for tempering substances which are placed on the heating plate in vessels. By means of a built-in stirring drive and with the help of a magnetic bar within the vessel, the

substances can be stirred at the same time. The mixing intensity is dependent on the motor speed and on the size of the magnetic bar. The RH basic KT/C and RH digital KT/C also has a contact thermometer connection in compliance with requirements of DIN 12 878 and an adjustable safety circuit.

Area of use

- › Indoor environments similar to that a laboratory of research, teaching, trade or industry area.
- › The safety of the user cannot be guaranteed:
 - if the device is operated with accessories that are not supplied or recommended by the manufacturer,
 - if the device is operated improperly or contrary to the manufacture's specifications,
 - if the device or the printed circuit board are modified by third parties.

Unpacking

Unpacking

- › Unpack the device carefully. Any damage should immediately be reported to the carrier (mail, rail or freight forwarding company).

Scope of delivery

- › An RH basic 2 or RH basic KT/C or RH digital KT/C
- › Operating instructions.

Commissioning

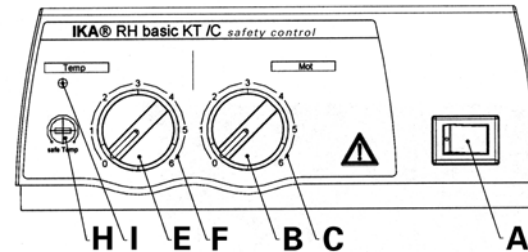
Check whether the voltage specified on the type plate matches the mains voltage available. The power socket used must be earthed (protective earth conductor contact). Before tuning on the machine, adjust the turn dial for controlling the speed so that it is against the left stop. If these conditions are met, the device is ready to operate after plugging in the mains plug. If these procedures are not followed, safe operation cannot be guaranteed and/or the equipment may be damaged.

Observe the ambient conditions (temperature, humidity, etc.) listed under Technical Data.


Starting

The heating and stirring features can be turned on and off by pressing the unit switch (A).

Version RH basic 2 and RH basic KT/C: The ready for operation and turned on states of the device are indicated by the green lamp in the unit switch lighting up.

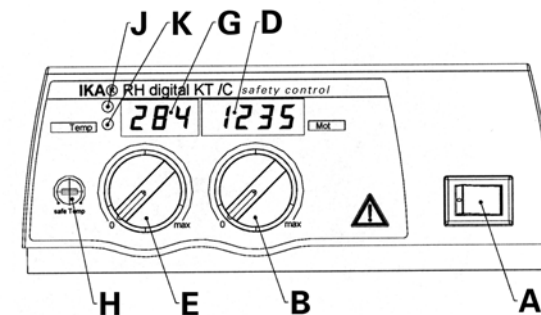Version RH digital KT/C: When you switch on the stirrer it will run a self-test and all the LED segments of the displays will light up for a few seconds. After this, the LED's (D; G) light up to indicate that the unit is powered up and ready for operation.

Function Stirring

The stirring rod in the glass beaker with a maximum length of 60 mm is driven by a permanent magnet located directly under the placing and heating surface. It allows magnetic field lines to pass through it. The permanent magnet is fastened directly onto the drive shaft of the motor.

The actual speed depends on the load and voltage. Note that fluctuations in the in the mains power voltage within the allowable tolerance and changes in viscosity of the materials to be stirred resulting from processes they undergo may also cause minor fluctuations in the speed.

Version RH basic 2 and RH basic KT/C: The speed of the motor can be adjusted

on the "Mot" dial (B) with a scale from 0 to 6 (C). The scale values from 0 to 6 on this dial correspond roughly to speeds ranging from 0 to 2000 rpm. The maximum speed of 2000 rpm is reached when the dial switch is turned all the way to the right and there is no load on the motor.

Version RH digital KT/C: Set the motor speed using the "Mot" knob (B); the knob has a graduated scale from 0 to maximum rpm. The motor speed is shown on the LED display (D). When the motor is stopped the speed "0" will be displayed. When the motor is operating, the display will show the current speed in the range 100 to 2000 rpm.

Function Heating

The temperature of the heating plate of the instrument is kept constant by a control circuit. In addition, it is monitored by a safety circuit. The required temperature sensor (a thermo-couple) is installed in the heating plate. The RH basic unit has a stainless steel surface plate with a heating output of 400 Watt. RH basic KT/C and RH digital KT/C units have a enamel-coated stainless steel surface plate with a heating output of 500 Watt.

Version RH basic and RH basic KT/C: The temperature of the heating plate can be adjusted on the „Temp“ turn dial (E) with a corresponding scale from 0 to 6 (F). The scale values from 0 to 6 correspond approximately to area temperatures up to a max. of 320 °C. When the setting is at 0, the device does not heat the medium up at normal room temperature. During the heating phase, the signal light (I) lights up orange (orange = power is being supplied to the heating plate; signal light off = no power being supplied)

Version RH digital KT/C: The temperature of the heating plate can be adjusted on the „Temp“ turn dial. When you change the knob setting, the LED display (G) will show the SETPOINT temperature and the warning light (K) will light up. After approx. 3 seconds the LED display will switch back to showing the ACTUAL

temperature and the warning light (K) will switch off. The temperature indication now constantly changes between actual temperature (25 sec) and target temperature (5 sec). When you select the SETPOINT temperature and as the unit heats up, the warning light (K) will remain off and the LED display (G) will show the ACTUAL temperature of the surface plate in the range from 0°C to +320°C. The red warning light (I) will light up while heating is in progress. (Red warning light ON = surface plate heating power supply is ON; red warning light OFF = surface plate heating power is OFF).

PLEASE NOTE!
The heating plate of the version RH basic KT/C and RH digital KT/C only heats up, if the contact plug has been plugged (the contact plug is undetachably fitted to the instrument), or if an ETC 1 and/or a contact thermometer has been connected.

Limiter of safety temperature

The safety temperature limiter makes it possible to adjust the maximum temperature of the heating plate (safety temperature). This ensures that the temperature of heating plate does not exceed the flame point of the material. The value to be set must be determined by trial and error and should generally fall about 20 K below the flame point of the medium. RH basic 2 devices have a safety circuit with a fixed setting that limits the temperature of a heating plate to 400 °C. RH basic KT/C and RH digital KT/C devices have an adjustable safety circuit. The maximum temperature of the heating plate can be adjusted within a range from 100 °C to max. 400 °C.

Adjusting the safety temperature limit "Safe Temp" (Vers. KT)

Version RH basic KT/C and RH digital KT/C: The "Safe Temp" set screw (H) is to the left of the "Temp" knob (E) on the front operator panel.

You can use the „Safe Temp“ set screw (H) of the safety temperature limiter to adjust the desired safety temperature. If the desired safety temperature is exceeded by a higher adjusted target temperature, or if there is a defect in the instrument controller, the heating turns off permanently in this state and **the stirring function continues to run at the speed that was set prior to the malfunction.**

- Using a straight-edge screwdriver, turn the "Safe Temp" adjustment screw (H) as far as it will go clockwise (to the right).
- Do not turn the screw beyond the natural right or left limit of rotation, as this will destroy the potentiometer.
- Use the "Temp" turn dial (E) to adjust the set-point temperature to desired safety temperature ("Safe Temp").
- Turn the "Safe Temp" set screw (H) slowly to the left until the heating function switches off and the signal light (I or J) goes off. Wait for about one more minute to ensure that the signal light (F) going off does not merely indicate the end of a switching cycle of the heater. If the signal light (I or J) turns on again within this time, you must repeat the procedure, slowly turning the "Safe Temp" (H) adjustment screw to the left again. When the signal light (I or J) does not light up again within one minute, the safety temperature is correctly set.
- After this, turn set "Safe Temp" adjustment screw (H) back slightly to the right (clockwise).
- After you have turned the unit switch (A) off and back on again, the device is ready for operation.

NOTE: When the "Safe Temp" (H) adjustment screw is set to the minimum level (as far as it will go to the left) the heating function is turned off.

Triggering of the safety circuit is shown in the following ways:
On the [RH basic 2](#) and [RH basic KT/C versions](#): the warning light (I) will flash two times per second. (Note: regular flashing indicates that the safety circuit has been triggered; this warning light remains permanently lit to indicate that a standard heating cycle is in progress).
On the [RH digital KT/C](#) version: the letter "-S-" will be shown on the LED display (G).

Controlling the Medium Temperature via ETC 1 or Contact Thermometer

The control of the medium temperature via ETC 1 or contact-thermometer should be preferred. Its advantages are: just a short heating-up phase after setting the rated temperature, practically no temperature drift and only a minor temperature waviness.

The rear of the instrument houses the diode bushing for connection the ETS-D or contact thermometer, or for plugging the contact plug.

The electronics of the instrument supply a test current that has to flow over plug pins 3 and 5 of the diode bushing in order for the heating plate to heat.

Safety contact thermometers

acc. to DIN 12 878 class 2 or acc. Gerstel are connected with a three-wire cable, the test current flows through the contact thermometer.

Safety function

If the test current is interrupted because of e.g. breakage of contact thermometer or falling out of the cable plug, the heating cuts off.

Contact thermometer without safety circuit

acc. to DIN 12 878 class 0. The instrument only heats if the test current circuit is closed by an electrical connection of the plug pins 3 and 5.

2-wire connecting cables:

Connect plug pins 3 and 5 of the instrument plug.

3-wire connecting cables:

Here the test current circuit can also be produced in the terminal head of the contact thermometer (connect plug pins 2 and 3).

A 3-wire cable with the required bridge is available (accessories).

The rated temperature is set on the contact thermometer by means of the rotary magnetic head, whereby the upper edge of the traveling nut is the reference point on the rated value scale.

The required surface temperature of the heating plat is preselected on the temp-rotary knob of the instrument.

If the temp-rotary knob is set to the maximally adjustable temperature, the heating-up takes place as rapidly as possible, yet the medium temperature overshoots the rated value set on the contact thermometer.

If the temp-rotary knob is set to twice the rated value, the result is good compromise between rapid heating-up and overshooting.

If the temp-rotary knob is set precisely to the rated temperature, the medium does not reach the rated temperature owing to a drop in temperature between heating plate and medium.

The rated temperature is setting on the ETS-D.

If you are using an ETS-D4, we recommend using the ETS-D4 in On/Off mode (two-point control) at higher medium temperatures.

Error codes

Version RH digital KT/C:

The LED displays (C and D) will start to flash and display a code when a fault is present. The fault codes and troubleshooting procedure are as follows:

| Error code | LED | Cause | Solution |
|-------------|-----|---|---|
| E 2 | G | Inner temperature too high | Switch off the unit and allow it cool down. Check that the ventilation slots on the bottom are not blocked and that cooling air can circulate freely. |
| - S - | G | Safety circuit is triggered | Reduce the temperature setpoint or reset the safety circuit where necessary. |
| E 2 | D | No pulse at the reading fork (faulty reading fork, faulty motor or motor blocked) | Switch off the unit. Trace the external cause for blocking the stirring drive. If no external cause can be determined, clean the unit and return it for repair. |
| E 6 | G | Contact plug not plugged in | Plug in contact plug |
| | | Breakage and/or short-circuit of the contact thermometers | Replace contact thermometer |
| Other codes | | | Switch off the unit. Clean and return it for repair. |

Version RH basic 2 and RH basic KT/C:

The warning light (I) flashes to indicate that a fault is present. To remove a fault condition, proceed as follows:

- Switch off the unit and allow it to cool down.
- Trace and remove the cause of the fault.
- Switch on the unit.

If the fault persists, clean the unit and return it for repair.

Mounting - Support rod H 16 V

The H 16 V support rod can be screwed into the threading on the right behind the housing as follows:

- Remove the plastic cap and washer from the support rod and screw off the hex nut.
- Place the washer and hex nut on the beginning of the threading on the housing and hold them in place.
- Screw the support rod into the nut and threading of the housing and counter-tighten the hex nut.

Maintenance and Cleaning

The device is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

Cleaning

- › For cleaning disconnect the mains plug!
- › Use only cleaning agents which have been approved by IKA to clean IKA devices.
These are water (with tenside) and isopropanol.
- › Wear protective gloves during cleaning the devices.
- › Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.
- › Do not allow moisture to get into the device when cleaning.
- › Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.

Version RH basic KT/C and RH digital KT/C

Information for Care and Maintenance of the Heating Plate with Technical Enamel Coating

The technical enamel makes the heating plate easier to care for and more resistant to acids and bases. Because of it, however, the heating plate is also more susceptible to extreme fluctuations in temperature and the force of impact. This can result in cracks forming or the coating flaking off.

Make certain that the bottom of the placing vessel is even, clean and dry. The bottom of the placing vessel must not have any sharp grooves, sides or edges. We recommend most strongly that you clean the heating plate regularly.

Spare parts order

- › When ordering spare parts, please give:
 - device type.
 - serial number, see type plate.
 - position number and description of spare part, see www.ika.com.
 - software version.

Repair

- › Please only send devices in for repair that have been cleaned and are free of materials which might present health hazards.
- › For repair, please request the “**Safety Declaration (Decontamination Certificate)**” from IKA or use the download printout of it from IKA website at **www.ika.com**.
- › If your appliance requires repair, return it in its original packaging. Storage packaging is not sufficient when sending the device - also use appropriate transport packaging.

Accessories

RS 1 set of stirring bars **H16 V** support rod
RSE PTFE stirring bar remover
Recommended IKA-stirrers
TRIKA-stirring bars: PTFE, triangular, length 55 mm
IKAFLON-stirring bars: PTFE diameter 8 mm, lengths 25, 30, 40 and 50 mm

For more accessories see **www.ika.com**.

Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from unintended use or insufficient care and maintenance contrary to the instructions in this operating manual.

Technical data

| | | RH basic 2 | RH basic KT/C | RH digital KT/C |
|--|------------|------------------------|--|---|
| Design voltage | VAC | | 220 ... 240 ±10% | |
| or | VAC | | 115 ±10% | |
| or | VAC | | 100 ±10% | |
| Design frequency | Hz | | 50/60 | |
| Motor power consumption | W | | 15 | |
| Motor power output | W | | 2 | |
| Speed range | rpm | | 100 ... 2000, infinitely variable setting | |
| Adjustment of the speed control | - | | rotary knob on the front | |
| Display of the speed control | - | | scale 0 ... 6 | LED-display four digit (0 - 2000 rpm) |
| Motor | - | | shaped pole motor | |
| Max. stirring volume (water) | ltr | 10 | 15 | |
| Heating plate | - | stainless steel 1.4301 | stainless steel 1.4301 with enamel coating | |
| Heating plate dimensions | mm | Ø 125 | 130 x 130 | |
| Heating power | W | 400 (230 VAC) | 500 (230 VAC) | |
| | W | 400 (115 VAC) | 500 (115 VAC) | |
| | W | 360 (100 VAC) | 450 (100 VAC) | |
| Surface temperature heating plate | °C | | RT ... 320 | |
| Temperature setting | - | | rotary knob on the front | |
| Temperature display | - | | scale 0 - 6 | LED-display four digit (0 °C ... +320 °C) |
| Contact thermometer connection | - | - | DIN 12 878 | |
| Infinitely variable safe temperature | °C | - | 100 ... 400 | |
| Fixed adjustment safe temperature | °C | 400 | - | |
| Permissible ambient temperature | °C | | + 5 ... + 40 | |
| Permissible relative humidity | % | | 80 | |
| Permissible on time | % | | 100 | |
| Screened for radio interference | - | | under VDE 0871 | |
| Protection class according to DIN EN 60529 | - | | IP 21 | |
| Contamination level | - | | 2 | |
| Overvoltage category | - | | II | |
| Protection class | - | | 1 (protective earth) | |
| Fuse | A | | 10AT Id. No. 2755400 | |
| Operation at a terrestrial altitude | m | | max. 2000 above sea level | |
| Dimensions | mm | | 168 x 220 x 105 | |
| Weight | kg | | 2.4 | |

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited