2000005144

UTTD_102019

IKA ULTRA -TURRAX® Tube Drive

IKA-Melko, Germany
Reg. No. 004343

Betriebsanleitung Ursprungssprache	DE	4
Operating instructions	EN	10
Mode d'emploi	FR	16
Руководство пользователя	RU	22
使用说明	ZH	28
Indicaciones de seguridad	ES	33
Veiligheidsinstructies	NL	34
Norme di sicurezza	IT	35
Säkerhetsanvisningar	SV	36
Sikkerhedshenvisninger	DA	37
Sikkerhetsanvisninger	NO	38
Turvallisuusohjeet	FI	39
Instruções de serviço	PT	40
Wskazówki bezpieczeństwa	PL	41
Bezpečnostní pokyny	CS	42
Biztonsági utasítások	HU	43
Varnostna opozorila	SL	44
Bezpečnostné pokyny	SK	45
Ohutusjuhised	ET	46
Drošības norādes	LV	47
Nurodymai dėl saugumo	LT	48
Указания за безопасност	BG	49
Instrucțiuni de securitate	RO	50
Υποδείξεις ασφαλείας	EL	51

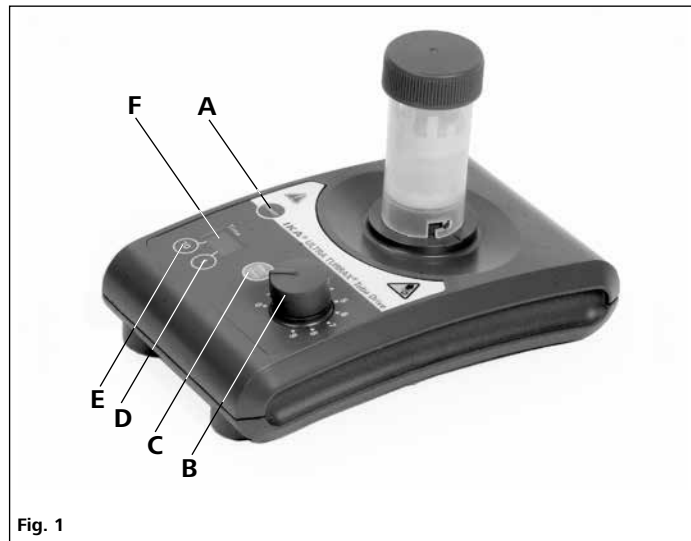
Fig. 1

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Contents

	Page
EU Declaration of conformity	10
Explication of warning symbols	10
Safety instructions	10
Unpacking	12
Correct use	12
Useful facts	12
Commissioning	12
Error message	14
Accessories	14
Maintenance	14
Warranty	14
Technical data	15

EU Declaration of conformity

We declare under our sole responsibility that this product corresponds to the directives 2014/35/EU, 2006/42/EC, 2014/30/EU and 2011/65/EU and conforms with the following standards or normative documents: EN 61010-1, EN 61010-2-051, EN 61326-1, EN 60529 and EN ISO 12100.

A copy of the complete EU Declaration of Conformity can be requested via: sales@ika.com.

Explanation of symbols

General hazard

This symbol identifies information **that is of vital importance for safeguarding your health and safety**. Disregarding this information can lead to health impairment and injuries.

This symbol identifies information **that is of importance for the technically correct functioning of the system**. Disregarding this information can result in damage to the device or to system components.

Risk of hand injury.

Risk of biohazard infection.

Safety instructions

For your protection

- **Read the operating instructions in full before starting up and follow the safety instructions.**
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- The device must be supervised at all times when in operation.

Note! Only suitable for diagnostic use in combination with **IKA IVD** Tubes (in according to IVD directive).

! DANGER

When handling infectious or toxic sample materials, do **NOT** touch the sample, mixture or waste with your hands. Wear gloves and lab coat and goggles if necessary.

! DANGER

The device must only be operated with a tube attached; with no tube there is a risk of injury from the rotating locating pin.

- Hearing protection must be worn when working with **IKA BMT-20** or **IKA BMT-50** tubes.

! DANGER

Wear your personal protective equipment in accordance with the hazard category of the medium to be processed. There is a risk of:

- splashing liquids
 - projectile parts
 - body parts, hair, clothing and jewellery getting caught.
- Set up the device in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
 - The feet of the device must be clean and undamaged.
 - The device is not suitable for manual operation.
 - The device may heat up when in use.
 - Check the device and accessories beforehand for damage each time when you use them. Do not use damaged components.
 - Ensure that the cover on the tube is screwed on tightly.
 - Ensure that the tube is firmly attached to the bayonet lock connector on the drive unit prior to operating the device.
 - The tube must only be attached and removed while the motor is stationary.

! WARNING

The **IKA** tubes must always be closed when the device is in operation. Switch off the device immediately if any material leaks from the tube. Clean the device.

- Always open the tube carefully after use as the media in the tube may heat up due to transfer of energy during operation, leading to pressurization of the container: risk of material spraying, protective equipment must be worn.
- The temperature of the material must not exceed 40 °C.
- Only use tubes approved by **IKA**.
- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- Do not use the device in explosive atmospheres, it is not EX-protected.
- With substances capable of forming an explosive mixture, appropriate safety measures must be applied, e.g. working under a fume hood.
- To avoid body injury and property damage, observe the relevant safety and accident prevention measures when processing hazardous materials.
- The device doesn't start up again automatically following a cut in the power supply.
- Safe operation is only guaranteed with the accessories described in the "**Accessories**" chapter.

For protection of the device

- The voltage stated on the name plate must correspond to the mains voltage.
- The device must only be operated with the original plug-in power supply unit.
- Protect the device and accessories from bumps and impacts.
- The device may only be opened by experts.

Unpacking

Unpack:

- Please unpack the device carefully.
- Any damage should be notified immediately to the shipping agent (post office, railway network or logistics company).

Delivery scope:

- **Tube Drive**
- Plug-in power supply with interchangeable adapters
- Removal hook
- Operating instructions
- A warranty card.

Intended use

Use:

	Application	Volume	Mode
▶ with Tube ST-20...	Mixing	2 - 15 ml	Seconds and minutes mode
▶ with Tube ST-50...	Stirring	15 - 50 ml	Seconds and minutes mode
▶ with Tube DT-20...	Dispensing	5 - 15 ml	Seconds and minutes mode
▶ with Tube DT-50...	Stirring	15 - 50 ml	Seconds and minutes mode
▶ with Tube BMT-20...	Crushing	2 - 15 ml	Seconds and minutes mode
▶ with Tube BMT-50...	Mixing wet/dry	15 - 50 ml	Seconds and minutes mode
▶ with Tube BMT-20-S-IVD ⓘ BMT-20-C-IVD ⓘ	Crushing Mixing wet/dry	2 - 15 ml	Seconds and minutes mode

ⓘ Note: for BMT-20-S-IVD/BMT-20-C-IVD:

- max. speed: scale 4
- max. Timer: 1 min.

Range of use:

Indoor environments similar to that a laboratory of research, teaching, trade or industry.

The safety of the user cannot be guaranteed:

- If the device is operated with accessories that are not supplied or recommended by the manufacturer
- If the device is operated improperly contrary to the manufacturer's specifications
- If the device is operated improperly contrary to the manufacturer's specifications.

Useful facts

Changes in viscosity and volume caused by processing of dispersion media may cause small fluctuations in the speed of the device.


Commissioning

Pay attention to the ambient conditions listed in the “**Technical Data**”.

Position the rotating knob (B) at the left stop before starting up.

If these conditions are met, the unit is ready for service when the mains plug has been plugged in.

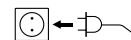
Switching on:

Press “Power” button (A).

Previous mode and time settings will be displayed e.g. :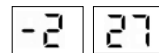


(Factory setting: -2)

Setting Timer (Changing of mode)

Switch off the device.

Minutes-mode

Press and hold "Start/STOP" button (C), and also press the "Power" button (A).

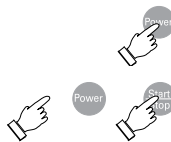Minutes-mode display:

The time that was previously set will be stored, e.g.:

Note: max. 29 min., speed range upto 4000 rpm.


Seconds-mode

Press and hold "Start/STOP" button (C), and also press the "Power" button (A).

Seconds-mode display:

The time that was previously set will be stored, e.g.: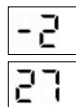

Note: max. 59 sec, speed range upto 6000 rpm.

Setting the time

The time that was previously set will be stored, e.g.:

(Factory setting: 00)

Set the tens digit using button (E).

Set the ones digit using button (D).

Attaching the tube

Fit the tube carefully and ensure it is properly attached.

Setting the speed

Set the desired speed in the range of 300–6000 rpm using the infinitely adjustable rotary knob (B).

Starting

Press "Start/Stop" button (C), while the preset countdown is running, timer display will blink.

An audible signal will be given when the countdown is complete.

Cancelling the countdown

Press button "Start/Stop" (C).

Switching off

Press button "Power" (A).

Error message

It is possible for the media to become lodged between the cutters of the DT-tubes. The unit will automatically change the direction of rotation up to three times in an attempt to free the cutters. The unit will switch off automatically if this is not possible.

If the device becomes blocked on three occasions within a certain period of time the following error message will occur:

⇒ Display:

⇒ An audible signal will be given.

Corrective action: Switch off the device, check contents, switch on the device.

Accessories

See more accessories on www.ika.com.

Maintenance

The device is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

→ For cleaning disconnect the mains plug.

Only use cleansing agents which have been recommended by **IKA**: Water (containing surfactant) and isopropyl alcohol.

- Wear protective gloves during cleaning the instruments.
- Electrical instruments may not be placed in the cleansing agent for the purpose of cleaning.
- Do not allow moisture to get into the instrument when cleaning.
- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with **IKA** that this method does not destroy the instrument.

Spare parts order:

When ordering spare parts, please give:

- machine type
- manufacturing number, see type plate
- item and designation of the spare part, see www.ika.com, spare parts diagram and spare parts list.

Repair

In case of repair the device has to be cleaned and free from any materials which may constitute a health hazard.

For repair, please request the "**Decontamination Certificate**" form **IKA** or use the download printout of it from **IKA** website: www.ika.com.

If you require servicing, return the appliance in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

Technical data

Power supply unit		
Input	VAC A Hz	100 ... 240 0.8 50 ... 60
Output	VDC Hz	24 24 W.LPS (limited power source) 50 ... 60
Protection class		2 (double insulated) 
Tube Drive		
Operating voltage	VDC mA	24 800
Power consumption, normal operation	W	20
Power consumption, standby operation	W	2
Power output	W	17
Drive		EC flat motor
Speed range:	rpm	300 ... 6000, infinitely adjustable
Mode 1		up to 4000
Mode 2		up to 6000
Speed setting		rotating/pressing knob
Speed display		scale 0 ... 9
Timer: Mode 1		1 ... 29 min
Mode 2		1 ... 59 sec
Display-timer		digital
Perm. ambient temperature	°C	+5 ... +40
Perm. rel. humidity	%	80
Protection to EN 60529		IP 20
Contamination level		2
Overvoltage category		II
Operation at a terrestrial altitude level	m	max. 2000
Dimensions (W x D x H) without tube	mm	100 x 160 x 40
Weight	kg	0.75

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IKA-Werke GmbH & Co.KG